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Table of Contents - Operating manual

1. Models
2. Technical specifications / Drawings
3. Intended application
4. General safety-related information / warranty
5. Transport
6. Setting up
7. Commissioning
8. Interior installations
9. Storage quantities / storage together
10. Ventilation
11. Marking
12. Maintenance / repairs
13. Operating instructions
14. Disposal

1. Models

- 1.1 CHS with wing doors
- 1.2 CHS with sliding doors
- 1.3 CHS with removable doors

2. Technical Specifications / Drawings

see Appendix 1

2.1 Design

Body: Powder coated sheet steel

Equipment: Height adjustable tray inserts or shelf inserts
Floor pan in accordance with § 19 WHG with Ü marking (Conformity Labelling)
according to the Steel trough guidelines (Stawa-R)

3. Intended application

storeLAB® chemical cabinets are used for storing water hazardous and / or poisonous substances in closed containers in workrooms as long as the storage is compatible with the protection of the employees.

4. General safety-related instructions / warranty

- The cabinets must be setup, operated and maintained in such a way that the safety of the employees is guaranteed against harmful vapours.
- The storage of combustible fluids, aggressive substances or materials that can lead to the development of fires and explosions due to spontaneously combustible or instable properties is forbidden.
- When operating the chemical cabinet, observe the respective national guidelines, directives and legislations that are used for this purpose.
- A functional check must be carried out before commissioning for the purpose of detecting damage caused by transport, shipment and incorrect installation!
- Guarantee / warranty exists according to our GTC. See also point 11.
- Please observe your obligation to exercise diligence - Damage must be eliminated immediately and leaks must be sealed!
- The opening range of the doors must be kept free and clean.
- Doors and drawers should only be kept open for supplying and removal - Keep the doors / drawers closed!
- Please observe the maximum storage quantity, loading of the shelf inserts, etc. according to Appendix 1 or according to the labels on the door / drawer.
- The chemical cabinets are not intended to be used for storing aggressive materials (e.g. acids and alkalis, corrosive gases) - otherwise there is a risk of corrosion to the interior equipment and the exhaust gas ! Use our special acids and alkalis cabinet for this purpose.
- Check the resistance of the cabinet surface before storing the storage goods (e.g. chemicals)
- The collecting equipment (floor pan) must collect 10% of the total amount stored above it but, at least the content of the largest single container. In water protection areas, 100 % of the storage quantity must be collected as long as the storage is permissible. (Applicable for Germany)
- The respective national guidelines must be fulfilled
- Suitable personal protective equipment must be worn when using substances or toxins harmful to health!

5. Transport

The storeLAB® chemical cabinets must only be picked up from the bottom side using suitable transporting equipment and must be transported vertically. Tilting when picking up and transporting must be avoided in all cases.

Always pick up the cabinet from the side or from the rear. Never pick up from the door side.

The pick up area of the barrow, forklift or similar must have at least the same floor space of the safety cabinet.

6. Setting up

The ground must be level and be able to bear the weight of the storeLAB® chemical cabinet when fully loaded.

The storeLAB® chemical cabinet must be setup horizontally and any safety equipment must be installed (e.g. wall holder with CHS-APO). This is the only way to ensure that the cabinet does not fall over.

The position of the actual location of use / operation must ensure that the chemical cabinet:

- cannot be damaged by vehicles,
- is not exposed to any sources of direct or indirect heat,
- is protected against moisture.

Suitable personal protective equipment must be worn when using substances or toxins harmful to health!

The location of the exhaust connections can be obtained in Appendix 1 (see sketches)

7. Commissioning

The chemical cabinet must be setup, operated and maintained in such a way that the safety of the employees is guaranteed.

With CHS-APO see assembling instructions (inside the cabinet door)

Suitable personal protective equipment must be worn when using substances or toxins harmful to health!

8. Interior installations

8.1 Tray insert

The supplied tray insert as well as the floor support are installed in the cabinet ex-factory.

For positioning the tray insert, lift the tray insert and pull horizontally from the cabinet. Pull the floor supports out of the bottom holes of the positioning strip and turn 90° upwards. Now pull the floor support standing horizontally from the top holes of the positioning strip. The assembly in the new position is carried out in the reverse order of the procedure previously described.

8.2 Extension tray

The supplied extension tray as well as the extensions with tray supports are installed in the cabinet ex-factory. For positioning the extension trays, pull these forwards, release the lock on the extensions and remove horizontally from the cabinet. Pull the extension trays with extension from the holes of the positioning strip and turn 90° upwards. Now pull the extension tray standing horizontally from the holes of the positioning strip. The assembly in the new position is carried out in the reverse order of the procedure previously described.

8.3 Mounting shelf

Already mounted in the cabinet

Observe the maximum loading capacity (see Appendix 1)!

If the shelf inserts are not mounted according to the assembling instructions, the maximum loading capacity cannot be ensured.

9. Storage quantities / storage together

The applicable national guidelines must be observed.

10. Ventilation

All chemical cabinets are provided with an exhaust air connection.. The supply air is drawn from the room. Please observe the national applicable guidelines, directives and legislations with regard to the prescribed air change and the exhaust air routing.

Operating storeLAB® chemical cabinets can also be carried out without technical exhaust air (in Germany, observe the TRGS here)

The respective national guidelines, directives and legislations must be observed in the process. The location of the exhaust connections can be obtained in Appendix 1 (see sketches)

11. Marking

Marking according to TRGS / ISO 3864/GHS-VO.

12. Maintenance / repairs

Daily operating inspections

Escaping liquids must be removed immediately and disposed off in a correct manner. Check the cabinet for mechanical damage, soiling and corrosion on a regular basis.

13. Operating instructions

When considering the operating manual and the operating conditions, the operator must create an operating instruction in a manner that can be understood and in the language of the employees operating.

The operator must make the operating instructions available at a suitable location at the place of work.

14. Disposal

The storeLab® chemical cabinet can be dismantled completely. The individual components can be dismantled and recycled. The national and local disposal guidelines must be observed. Parts of the chemical cabinet or the entire cabinet must not be disposed of as bulk rubbish or domestic waste for the purpose of preserving resources.

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to which this declaration relates is in conformity with the relevant provisions of the following standards or other
normative documents.

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**EG-Maschinenrichtlinie 2006 /42/EG, L'157/65
vom 9.6.2006**
EEC-Machinery Directive 2006/42/EEC

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Applied national standards (germany) DIN 4844-1 Safety colours and safety signs

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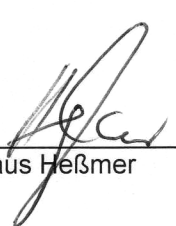
The operation company is responsible for installing the products in accordance to national regulations which may apply in respect to the installation, use and also for regular inspections.

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LA CONT Umwelttechnik GmbH